

Method Sheet 12

Plating macrophages for cytokine production assay

Overview

This method sheet explains how to seed 96-well plates with adherent macrophage cell-lines, such as RAW, J774 or THP-1, for the purpose of screening plant extracts for anti-inflammatory molecule discovery.

Reagents

- A growing culture of a mammalian monocytic or macrophage-like cell-line (e.g. RAW 264.7, J774, J774A.1, THP-1)
- Optional: If using THP-1 cells, phorbol 12-myristate 13-acetate (PMA) solution at 1 mM in ethanol or DMSO
- RPMI with L-glutamine, 10% FCS and antibiotics

Equipment

- Sterile, clear plastic 96-well microplate(s) with lids
- Multichannel pipette (8- or 12-channel) capable of dispensing 100 - 200 μ l
- Sterile pipette tips (compatible with the multichannel pipette)
- Sterile reservoir(s) for cell suspensions to set up the plate
- Tissue culture suite with a Class 2 biosafety cabinet and an incubator capable of maintaining a 5% CO₂ atmosphere at 37°C
- Waste container for used tips

Plating cells (RAW or J774 macrophages)

- 1) One day before challenge, suspend adherent RAW or J774 cells from a T75 flask using a sterile plastic scraper (trypsin / EDTA is not necessary for these cell-lines).
- 2) Pellet the cells by centrifugation (e.g. 300 g for 5 minutes).
- 3) Discard the supernatant and resuspend the cell pellet in 10 ml RPMI / 10% FCS.
- 4) Count the viable cells using a haemocytometer or automated cell counter.
- 5) Adjust the cell suspension to a concentration of 4×10^5 cells per ml in a total volume of 12 ml (enough for one 96-well plate, multiply if you intend to prepare more plates).
- 6) Pour the cell suspension into a suitable sterile plastic reservoir.
- 7) Use a multi-channel pipette to aliquot 99 μ l of cell suspension into every well of a 96-well plate.
- 8) Incubate the cells at 37°C in a 5% CO₂ atmosphere overnight (12-24 hours) before challenge.
- 9) This volumes given in this method are sufficient to prepare one 96-well plate, multiply appropriately if you intend to prepare more plates.

Plating cells (THP-1 macrophages)

- 1) Three days before challenge, suspend adherent RAW or J774 cells from a T75 flask using a sterile plastic scraper (trypsin / EDTA is not necessary for these cell-lines).
- 2) Pellet the cells by centrifugation (e.g. 300 g for 5 minutes).
- 3) Discard the supernatant and resuspend the cell pellet in 10 ml RPMI / 10% FCS.
- 4) Count the viable cells using a haemocytometer or automated cell counter.
- 5) Adjust the cell suspension to a concentration of 4×10^5 cells per ml in a total volume of 12 ml.
- 6) Pipette 199 μ l of RPMI / 10% FCS into a separate, sterile 1.5 ml microtube and supplement with 1 μ l of the 1 mM PMA stock to yield a 5 μ M sub-stock.
- 7) Supplement the 12 ml of THP-1 cells with 120 μ l of the 5 μ M PMA solution and mix well (the final concentration of PMA in the cell suspension is now 50 nM).
- 8) Pour the cell suspension into a suitable sterile plastic reservoir.
- 9) Use a multi-channel pipette to aliquot 99 μ l of cell suspension into every well of a 96-well plate.
- 10) Incubate the cells at 37°C in a 5% CO₂ atmosphere overnight for 3 days before challenge.
- 11) These volumes given in this method are sufficient to prepare one 96-well plate, multiply appropriately if you intend to prepare more plates.

Challenging cells with *Phytotitre* extracts and LPS

- 1) Perform this step 12-24 hours after plating J774 or RAW macrophages, or 3 days after plating THP-1 macrophages (follow Method Sheet 13).

Typical timetables for plating, challenging and harvesting tasks

Day	Task (J774 or RAW cells)	Task (THP-1 cells)
Mon	Split / passage stock culture and prepare 96-well plate(s)	Split / passage stock culture and prepare 96-well plate(s) with PMA
Tue	Challenge cells in 96-well plate(s) with extracts and LPS, collect and freeze supernatants 3 hours later	-
Wed	-	-
Thu	-	Challenge cells in 96-well plate(s) with extracts and LPS, collect and freeze supernatants 3 hours later
Fri	Split / passage stock culture	Split / passage stock culture

Notes

- The PMA step is necessary to promote differentiation of the THP-1 pro-monocytic cell-line from its typical suspension form into an adherent, macrophage-like form that is also more responsive to LPS.
- Remember to supplement the stock bottle of RPMI with L-glutamine if this is not present in your variant of RPMI medium.
- Never add PMA to the stock flask of growing THP-1 cells, this would halt their proliferation permanently.
- PMA treatment is not necessary for the J774 or RAW macrophage cell-lines as they are already adherent and sensitive to LPS treatment.
- PMA is barely soluble in water, so ethanol or DMSO is used as the solvent for the high concentration stock solution.
- It is possible to perform more than one experiment per week with the THP-1 cells, by plating on Friday and challenging on Monday.

Disclaimer: These method sheets and other resources are provided for educational purposes only. The user's University Supervisor remains the Principal Investigator and the sole party responsible for the safe conduct, risk assessment, and ethical oversight of all laboratory work. Caithness Biotechnologies Ltd. accepts no liability for any injury, loss, or damage resulting from the application of the advice or protocols provided herein. Copyright © 2026, Caithness Biotechnologies Ltd. All Rights Reserved.