

Method Sheet 14

Measuring TNF- α secretion using the L929 cell bioassay

Overview

This method sheet explains how to perform the L929 cell bioassay to quantify the pro-inflammatory cytokine tumour necrosis-factor (TNF)- α , with the aim of screening a natural product library for anti-inflammatory molecule discovery. L929 cells are a mouse fibroblast cell line that become sensitised to the cytotoxic effects of TNF- α when exposed to actinomycin-D. Sensitised L929 cells are challenged in 96-well plates with supernatants from macrophages previously stimulated with bacterial endotoxin (LPS) to release pro-inflammatory cytokines. In this setting, TNF- α present in the macrophage supernatant causes L929 cell death, which is measured by crystal violet dye staining of remaining adherent, living L929 cells. The concentration of TNF- α in the tested supernatant is inversely related to absorbance measured at 540 - 600 nm. Natural extracts that inhibit the loss of crystal violet staining in this assay may therefore contain anti-inflammatory molecules. This method, based on the technique of Delahooke *et al.* (1995)¹, can be a more cost-effective alternative to ELISA for quantitation of TNF- α production, particularly in multi-plate screening programmes.

Reagents

- A growing culture of the L929 mouse fibroblast cell-line
- Stock actinomycin-D (Δ toxic, handle with care) at 1 mg/ml in DMSO or ethanol (store at -20°C)
- DMEM / 10% FCS (5% FCS works equally well at all stages of the assay)
- Crystal violet reagent (0.2% crystal violet in 20% methanol, see Notes section for how to prepare)
- 20% acetic acid in dH₂O (see Notes section for how to prepare)
- Optional if standard curve desired for modified experiments: 10,000 IU/ml TNF- α in PBS (store in small aliquots at -20°C)

Equipment

- Sterile, clear plastic 96-well microplate(s) with lids
- Sterile pipette tips (compatible with each pipette)
- Sterile reservoir(s) to dispense sensitised L929 cells
- Multichannel pipette (8- or 12-channel) capable of dispensing 1 - 10 μ l
- Multichannel pipette (8- or 12-channel) capable of dispensing 100 - 200 μ l
- Tissue culture suite with a Class 2 biosafety cabinet and an incubator capable of maintaining a 5% CO₂ atmosphere at 37°C
- Squeezable wash bottle dispensing distilled water (H₂O)
- A large plastic tray to collect waste runoff from plate washing
- Waste stream for halogenated chemicals

- Microplate reader capable of measuring absorbance of 96-well plates at 570 nm (or alternatively, any wavelength between 540 nm and 600 nm)

Day 1 (morning) - Seeding plate(s) with L929 cells

- 1) Stock L929 cells are cultured in DMEM supplemented with 10% FCS and penicillin / streptomycin.
- 2) At any time in the morning, split / passage L929 cells from the stock flask by adding trypsin / EDTA and incubating for 2 - 5 minutes at 37°C.
- 3) Note that cell clumping limits the performance of this assay, so it is helpful to ensure that trypsinisation proceeds for long enough to yield a suspension of single cells at this stage.
- 4) Pellet the cells by centrifugation (e.g. 300 g for 5 minutes).
- 5) Discard the supernatant and resuspend the cell pellet in 10 ml DMEM / 10%FCS.
- 6) If cell clumps remain visible by microscopy, pipetting up and down gently several times can help assist the process of disaggregating towards a suspension of single cells.
- 7) Use ~1 ml to seed a new stock flask, then count the remaining cells using a haemocytometer or an automated cell counter.
- 8) Prepare 12 ml of L929 cells at a concentration of 4×10^5 cells/ml in DMEM / 10%FCS (this is sufficient for one 96-well plate, multiply accordingly if preparing more than one plate).
- 9) Supplement the L929 cell suspension with 1:1,000 volume of 1 mg/ml stock actinomycin-D solution (e.g. add 12 μ l actinomycin-D to 12 ml of cell suspension).
- 10) ⚠ Note that actinomycin-D is very toxic - it must be handled with care, ideally with double gloves.
- 11) Place the cap on the tube and mix well by gentle inversion several times.
- 12) Pour the L929 cell suspension into a sterile plastic reservoir.
- 13) Use a multichannel pipette to transfer 90 μ l of this suspension into every well of a 96 well tissue culture plate.
- 14) Incubate the cells at 37°C for between 3 and 6 hours to allow the cells to adhere.

Day 1 (afternoon) - Challenging plate(s) with macrophage supernatants

- 1) At least 3 - 6 hours after plating the L929 cells, defrost the plate containing macrophage supernatants (prepared as per Method Sheet 13) in a 37°C incubator (takes about 20 minutes to thaw completely).
- 2) Remove the plate containing the L929 cells from the incubator to a Class 2 biosafety cabinet.
- 3) Align the L929 cell plate next to the macrophage supernatant plate, with well A1 on both plates at the top left position.
- 4) Use a multichannel pipette to carefully decant 10 μ l of supernatant from each well of the macrophage supernatant plate, and place the liquid directly into the respective wells of the plate containing L929 cells.


- 5) It is vital that the monolayer of cells at the bottom of the well is not scratched by the pipette tip during this procedure, so do not push the tips too far below the surface of the liquid.
- 6) Replace tips with fresh ones in between each column of supernatants transferred.
- 7) Return the plate containing L929 cells to the 37°C incubator and culture overnight.

Day 2 - Staining L929 cells with crystal violet dye

- 1) Cell culture methods preceding the staining step should be performed in a sterile environment with attention to sterile technique.
- 2) However, after the cells have been challenged, the staining part of the crystal violet assay does not need to be performed in a sterile environment - the following steps can be performed on an open bench in any laboratory with a suitable chemical waste stream.
- 3) Note that this assay can be messy, so it may be helpful to avoid stains from accidental spills by covering the bench surface with protective paper (such as Benchkote).
- 4) It is essential to wear appropriate PPE for this protocol, including lab coat, gloves and safety glasses.
- 5) Aliquot 6 ml crystal violet solution (see notes) per microplate to a suitable reservoir.
- 6) Remove all medium from the wells by inverting vigorously over the plastic tray, then patting dry with paper tissue.
- 7) Adjust a multichannel pipette to dispense 50 µl.
- 8) Firmly press the pipette into a row of sterile pipette tips (8 or 12, depending on the type of pipette), ensuring all are securely attached.
- 9) If the tips are loose or fall off during use, use gloved finger and thumb to pull up and seat firmly each tip individually, being careful to not get dye on your glove.
- 10) Push the plunger on the pipette down to the first stop (not all the way to the second stop).
- 11) Insert the tips into the liquid in the reservoir, ensuring all tips are below the surface of the liquid.
- 12) Slowly release your thumb to allow the plunger to return to the top position.
- 13) Look carefully across all the tips to ensure the level of liquid is the same in each tip, if not, dispense the suspension back into the reservoir and try again.
- 14) Likewise, if there are any large air bubbles in any of the tips, dispense back into the reservoir and try again.
- 15) Move the pipette to the plate containing cells and dispense all the liquid into the wells of the next available column, pipetting past the first stop all the way to the second stop of the plunger this time.
- 16) Repeat this process until every well in the plate has received crystal violet reagent.
- 17) Replace the lid on the plate and move the plate gently backwards and forwards to ensure the reagent covers completely the bottom of every well.
- 18) Incubate the plate at room temperature for **15 minutes**.
- 19) Gently wash all crystal violet solution from every well using distilled water from a squeezable wash bottle (tap water is also acceptable for this step).

- 20) Invert the plate over a large plastic waste tray until all liquid has been expelled.
- 21) Pat the plate dry after each wash by inverting the plate onto tissue.
- 22) Repeat the process of filling the wells gently with water, inverting the plate to remove the water and then drying with paper tissue, three more times.
- 23) The plate is correctly washed when little or no more blue reagent appears on the tissue.
- 24) Use a multichannel pipette to add 100 µl of 20% acetic acid solution to every well.
- 25) Pat the plate gently at the side to solubilise the cells until the colour in each well reaches visible homogeneity (about 30 seconds).
- 26) Use a microplate reader to measure absorbance of every well at 570 nm (filters between 550 and 600 nm will also work well for this purpose).
- 27) Retrieve the results from the microplate reader software for later analysis.
- 28) Empty the liquid contents of the plate into the waste tray, and discard the waste plastic to a suitable dry waste stream.
- 29) Discard the contents of the waste tray into a waste bottle that has been designated for the disposal of chlorinated chemical waste (often found in a fume cabinet).
- 30) Wipe the waste tray dry with tissue for re-use.
- 31) Clean up any spills by first wiping up any excess with tissue, then spraying the stain with 20% acetic acid, wiping with tissue, then spraying with 100% H₂O, then wiping again, repeat until all stain is removed.
- 32) Proceed to analysis of the crystal violet staining data to identify inhibitors of LPS-induced TNF-α production (follow Method Sheet 21).

Notes

-  Actinomycin-D is very toxic - it must be handled with care, ideally with double gloves.
- It is essential to wear appropriate PPE for this protocol, including lab coat, gloves and safety glasses.
- Be very careful not to get this stain on the bench, your skin or clothes - the stain can be messy and is very difficult to remove.
- If you do spill crystal violet stain on a surface, wipe up the excess with tissue, then spray with 20% acetic acid, wipe up with tissue, then with 100% H₂O, wipe again with tissue, repeat until all stain is removed.
- The methanol in the crystal violet reagent is the fixative that hardens the adherent cells.
- Methanol is toxic, volatile and flammable - prepare the working solution in a fume cupboard.
- Dead cells are washed away during the washing steps as they are non-adherent.
- The acetic acid in the resuspension reagent solubilises the crystal violet dye, allowing a uniform colour to be measured by a microplate spectrophotometer.
- Be careful to not wash the cells too vigorously - point the nozzle of the wash bottle at an angle away from the bottom of the wells by holding the plate at a 45° angle from vertical, and squeeze gently.

- Excessive pressure during the washing step will result in loss of adherent cells and poor reproducibility.
- It is possible to modify this assay to quantitate the bioactivity of TNF- α in macrophage supernatants relative to a recombinant TNF- α control, see Method Sheet 43.
- ¹ Method reference: Delahooke DM, Barclay GR, Poxton IR. Tumor necrosis factor induction by an aqueous phenol-extracted lipopolysaccharide complex from *Bacteroides* species. *Infect Immun* 63:840-6 (1995)

Disclaimer: These method sheets and other resources are provided for educational purposes only. The user's University Supervisor remains the Principal Investigator and the sole party responsible for the safe conduct, risk assessment, and ethical oversight of all laboratory work. Caithness Biotechnologies Ltd. accepts no liability for any injury, loss, or damage resulting from the application of the advice or protocols provided herein. Copyright © 2026, Caithness Biotechnologies Ltd. All Rights Reserved.